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Safety Precautions for Using a Forced Vacuum Drying Oven

NO. 035

Prohibitions

1. Strictly prohibited to heat first and then evacuate/vacuumize, to avoid cracking of the observation window and damage to the vacuum pump.
2. It is strictly prohibited to place highly corrosive, flammable, or explosive reagents, as well as flammable items such as plastic baskets and paper, inside the oven.
3. It is strictly prohibited to heat non-pressure-bearing sealed containers.
4. Do not stack items around the drying oven as this may affect heat dissipation. Flammable or explosive chemicals are strictly prohibited in the vicinity.

Operating Guidelines

1. Vacuum Evacuation Procedure:
 - Ensure that the balance port and vacuum valve are closed.
 - Turn on the vacuum pump first, and then open the vacuum valve.
 - After the vacuum level reaches the desired value, be sure to close the vacuum valve first, and then turn off the pump, to prevent back flow of pump oil.
2. End of drying procedure: Allow the temperature to cool down to below 60°C, then slowly open the balance port to return to atmospheric pressure, and subsequently remove the sample.
3. For powder sample drying: The sample container shall provide adequate coverage to prevent powder from splashing or becoming airborne during evacuation, while maintaining sufficient ventilation openings to ensure proper air/gas flow within the chamber.
4. After use, be sure to cut off the power supply and promptly clean out any items and residual stains inside the chamber.
5. For overnight operation of the drying oven, prior approval for overnight experimentation is required.
6. When the equipment triggers an over-temperature alarm, the power must be cut off immediately. After troubleshooting the fault, the equipment can be restarted.